

We supply:

- Living materials
- Culture media
- Frozen materials
- Preserved materials
- Skeletons
- Microscopes
- Wall charts
- Stains and Indicators
- Seeds
- Books
- Chemistry Models
- Anatomical Models
- Earth Science
- Environmental
- Forensic Equipment

and lots more on our website

www.southernbiological.com

Algal balls

The algal balls will arrive stored in a container of distilled water.

The *Chorella* within the balls is alive and therefore needs light to stay active. We recommend that you leave the balls in the container that they arrive in, and that you place this vial in a well lit area until required. Do not place in direct sunlight, nor expose to hot light sources, as the heat will be detrimental.

It will also help to loosen the lid of the container to allow air access. However, take care to protect the container from being knocked and subject to leaking.

For best results, use the algal balls as soon as possible, but in any case within 2 weeks of receipt.

Catalogue Code:
L1.60 and L1.61

***Supporting Products for
Chlorella cultures:***

L1.60 Algal photosynthesis kit

SI20 Hydrogen Carbonate Indicator

SI21 Hydrogen Carbonate, single set of standard solutions. (pH7.6 - 9.2 in increments of 0.2 pH units)

SI22 Hydrogen Carbonate, volume set of standard pH solutions, vials and Concentrate Hydrogen Carbonate indicator. (pH7.6 - 9.2 in increments of 0.2 pH units)

Ref: L1.60, L1.61 19/02/2016

1/44 Rushdale Street, Knoxfield VIC 3180 or PO Box 9227, Scoresby, VIC 3179, Australia
ph: 03 9753 3896 fax: 03 9753 3959

Email: sales@southernbiological.com Web: www.southernbiological.com
A Division of Cogitamus Pty Ltd ACN 097 468 954 ABN 25 838 926 933